

ticl



## Deep cooled scientific SWIR camera

# SIRIS

#### InGaAs sensor

> FPA 640x512, 0.9- 1.7 μm

#### Ultra-low Read-Out Noise

> below 5e-

#### Vibration and Cryogenic-Free Cooler

> selectable T° down to 77 K

#### Ultra-high Dynamic Range

> in Lin/log mode : > 140 dB

#### High speed sensor


> up to 200 fps full frame

## SIRIS

#### High-performance SWIR Camera

Short-wave InfraRed Imaging System

SIRIS is the most versatile and best performing scientific SWIR InGaAs camera on the market, combining ultra low noise, ultra large dynamic range and high speed. The camera is field proven on many observation missions at the T1M telescope of the Pic du Midi over the past couple of years. Only the most reliable components are integrated, like a vibration-less (below 1um) 200.000hrs MTBF Stirling cooler or a high-grade FPGA embedded platform for the most demanding applications you can imagine.

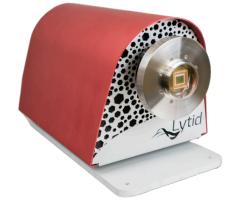


With SIRIS, SWIR photons have no where to hide from your eyes!

### Noise reduction with Non Destructive Read Out

With SIRIS, the longer you expose, the lower the noise! Using Non Destructive Read Out (NDRO) you can read the accumulated charges during an exposure without destroying them, and therefore achieve efficient readout noise reduction. SIRIS's exclusive NRDO read-out noise reduction allows to drastically reduce the read-out noise to *less than 5e-* using intra-exposure frame averaging. This unique feature on the market combined with a deep-cooled sensor (77-150K) can let you acquire shot-noise limited images in ultra-low light applications.

**M15** : with **100 NDROs** during a 1s exposure (right). The measured noise is 14e-, resulting in a reduction of approximately **6 times** when compared with standard exposure (left). This is equivalent of **2 magnitude** detection gain between the two images.

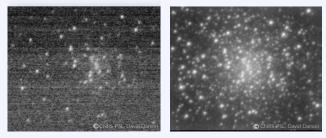

STANDARD

NDRO



SIRIS



HUBBLE SPACE TELESCOPE IN VISIBLE AND IR

A close view of a filed in the very center of the **ORION NEBULAE** (with the trapeze). Stack of exposures of 2s from our SIRIS SWIR Camera (left) with NDRO noise reduction . On the right, the same field taken by the SWIR camera of the Hubble Space Telescope.

**URANUS** : Stack of 400ms subexposures with **NDRO** noise reduction. Satellites, pole clouds, and **rings** can be seen. Same image, inverted confirms rings observation.







ORIGINAL

INVERTED

Zoom on URANUS




## Dual read-out mode : -CTIA -Lin/log

SIRIS provides two read-out modes, full linear (CTIA) and linear/logarithmic for class-leading >140dB dynamic range. Three adjustable gain levels ensure flexibility to suit broad variety of illumination conditions. Long exposure time up to one hour is achievable, and selectable region-of-interest on the detector allows exceptional frame rate values from 200fps full frame up to more than 10k fps ROI.

**SATURN** in Lin/log : With just one frame of 4s exposure time, on the right the same image as the left one but with viewing levels streched and turn on negative to better see low signals objects.



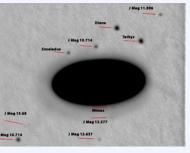

ORIGINAL

INVERTED



1175 nm



1275 nm


be observed by SIRIS with stack of 200ms exposure time at wavelengths of **1175nm** (50nm FHWM) filter and **1275nm** (50nm FWHM) filter with **no NDRO**.

**VENUS in CTIA** : Dark side of Venus could

### Highest dynamic range, lowest read noise ever achieved

Thanks to our new ultra-HDR (uHDR) method combine with the use of NDRO, the best ultra-high dynamics range images could be achieved. This is possible with the use of CTIA mode (ie full linear), on the same acquisition of very high signals and very low ones.


JUPITER with METIS and AMALTHEA or THEBE on the same raw images. The two images are from a stack of 32 1s subexposure with NDRO use for noise reduction combined with uHDR. The magnitude of Jupiter is -2.7 and Metis is 17.5. Magnitude difference of 20.3 on this image with only use of NDRO during 1s sub-exposure frame.

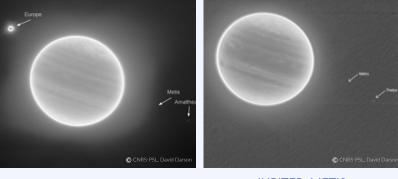

JUPITER, METIS, AMALTHEA JUPITER, METIS, THEBE

All pictures has been aquired with SIRIS on 1m Telescope at Pic du Midi



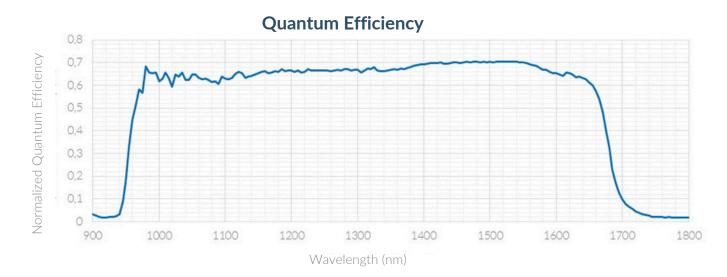

	ifications SIRIS
Detector	
Туре	InGaAs
Resoluti	on 640 x 512
Spectral	Response 0.9 – 1.7 µm
Pixel siz	e 15 μm
Dual-mo	ode sensor CTIA-linear Lin/Log
Performan	ce
Dark sig	nal <10e-/s @ 150 K
Gain	3 pixel-gain levels
Read-ou	t-modes Standard & NDRO
Read-ou	t-Noise <50e- lin mode, high gain <5e- NDRO
Well dep	th 300ke-, lin mod, low gain
Dynamic	s >140dB, lin/log
Digitizat	on 16-bit
Shutter	Global & NDRO
Region-c	f-interest ROI on detector, configurable
Frame-ra	200 fps full frame > 10 000 fps with ROI
Exposure	e-time From 1 μs to 1h
Trigger	Trig. In and Out (to 10ns)
Commur	nication Camera link
Cooling	300 K-77 K, cryocooler
Dimensio	ons 41 x 25 x 23 cm

Lytid SAS 10 rue A. Domon et L. Duquet 75013 Paris - FRANCE ☆ sales@lytid.com ⊕ www.lytid.com

US & Canada: Boston Electronics www.boselec.com thz@boselec.com +1.617.566.3821

